



# **13 THINGS TO KNOW ABOUT MOISTURE ANALYZERS**

*by Rachel Kohn, Ph.D.*

Tovatech LLC









---

**You Mentioned Display. What is Displayed?**

Backlit LCDs provide a wealth of information, the extent of which depends on the particular model. Examples include current weight, current moisture content in %, the drying mode (see below), the switch-off mode (see below), the previous drying time, the current temperature, and an indicator that the drying process is active. As suggested above data can be transferred via an RS232 interface to a printer or computer for record keeping. Units are available with software delivering a real-time graphic display of the moisture loss profile to your computer.



**How is Heat Applied to Samples?**

Halogen or infrared heaters are commonly used. These are mounted on precision balances. The instrument can be programmed to conduct the analysis in accordance with procedures recommended by industry trade associations and government agencies. As to the difference, halogen is much faster.

Temperature is generally maintained within 1°C of the selected setting. While most analyses involve increasing the temperature to a maximum of 160°C or less, in special cases higher temperatures may be necessary to achieve complete drying.

Note that there are other methods including microwave and the Karl Fischer method. Microwave drying instruments are somewhat larger and significantly more expensive than halogen-heated moisture analyzers. The Karl Fischer titration method, while accurate, uses hazardous solvents, requires skilled technicians and that the water in the sample be accessible to the reagent in order for the method to work. The method detects only water, whereas thermogravimetric analyzers detect any volatile substance in the sample.

## **What are Drying Modes or Profiles?**

Not all substances react equally to a moisture analyzer. Because of this some models offer a programmable heating profile. The most commonly used mode is standard drying, which is suitable for most substances.

The soft or gentle drying mode is recommended for samples that do not tolerate the full power of the halogen heater. This helps protect sensitive samples from decomposing or forming a protective “skin.” In the soft drying cycle samples are heated evenly from the outside to the inside by heat convection.

For samples with moisture content between 5% and 15% a fast drying cycle can be programmed into the unit. With this setting the halogen heater quickly raises the temperature to approximately 40% above the set temperature (called preheat) to rapidly drive off the bulk of the moisture and then returns it to the set temperature to approach the end point slowly

Step drying can be used in the same way as fast drying primarily for samples with more than 15% moisture. It can also be used for measuring two different types of moisture in the same sample, such as surface moisture and bound moisture. Individual increments of time and temperature are programmed into the moisture analyzer.



### **How Important is Sample Preparation?**

Very important in terms of sample surface area, size, weight and placement on the moisture analyzer sample pan. Here are some suggestions:

- Samples should be representative of the bulk material being tested.
- Samples should be tested immediately after being taken from the bulk material. Otherwise they could give up to or absorb moisture from the local environment.
- Small particles will analyze quicker because of increased surface area. Grains and beans, for example, should be crushed before testing.
- Samples should be spread thinly and evenly on the sample pan. Samples with high moisture or oily content – such as honey, milk, and body lotion – should be placed between fiberglass filters so they spread evenly across the sample pan. This not only increases surface area to speed the analysis but also prevents spattering inside the unit.

**What is the Correct Sample Size?**

The optimum sample weight is usually established after several trials. Moisture analyzer operating manuals may suggest weights for specific products but these are only suggestions. In general the sample should be sufficient to thinly cover the surface of the sample tray and evenly distributed to avoid lumps or peaks. Large sample sizes do not equate to more accurate readings except as noted below. The moisture balance operating manual will specify its maximum capacity.

Tip: Low moisture values such as below 1% are difficult to gauge, and require larger samples. As an example, for a 1% moisture content you may need a 2 gram or more sample, but for a sample with a 0.1% estimated moisture content you may need a 20 gram or more sample. Lower moisture samples require a more sensitive moisture analyzer. This is determined by its readout.

**What are Switch-Off Criteria?**

Switch-off criteria are selected by the operator depending on the samples under analysis and the required accuracy. Three methods of shutoff are common to most moisture analyzers: (1) automatically when drying is complete and the dry weight is stable, (2) after a specified time, or (3) manually. A stable dry weight is established when the weight loss per unit of time falls below a specified value.

**What is Calibration and Why is it Important?**

As with all precision balances accuracy is a paramount requirement for a moisture analyzer. Accuracy is set and maintained by initial calibration and periodic recalibration, the latter of which is performed based on schedules contained in a company's operations manual.

In the case of moisture analyzers, two calibration processes should be performed: one for weighing accuracy and one for temperature accuracy.

"External calibration" is the process used to check the weighing accuracy of most moisture analyzers. This is performed by using the recommended adjustment weight (such as 50 g) for the particular moisture balance being tested. Calibration procedures are outlined in the operating instructions accompanying various models of moisture analyzers.

Temperature calibration kits are available to check the accuracy of the heaters. Temperature tests are conducted at two points and take longer than the weight calibration exercise. Instructions are provided to perform the correcting adjustment(s) if there is a discrepancy between the temperature displayed on the probe and the analyzer.

## How are Moisture Balances Maintained?

Maintenance instructions are provided in the user manual. Here is a general summary:

First of all, position your moisture analyzer in a protected area away from drafts, dust, temperature extremes and vibration. Remember that every time you move it you must recalibrate it.

Use a soft brush or gentle vacuum to immediately remove samples that fall off the sample plate. Use extreme caution to ensure no liquid penetrates the device.

For general maintenance unplug the unit and remove accessory parts such as the sample dish, the dish retainer, the sample dish holder, the draft shield and its associated base plate. These can be individually washed using soft cloth dipped into water containing a mild liquid dishwashing soap and wrung out. Do not use solvents or other aggressive cleaning agents.

After cleaning wipe the components with a cloth dampened with fresh water then dry them with another soft cloth.

The same cautions apply to the body of the moisture balance.

The moisture balance temperature sensor is an extremely delicate component. Along with the halogen lamp, it is located in the lid. Cleaning is done in the same fashion as for the base of the unit. Avoid touching the halogen lamp. If you must touch the lamp use lint-free gloves to avoid transferring oil from your skin to the lamp surface.

**Do you have a Short List of Specifications to Check Out?**

Maybe not so short but here is a quick summary of things to look for when comparing moisture analyzers:


1. Readout (also called resolution or readability). This is the smallest unit that can be displayed on the panel and is presented in grams and percent moisture. Many moisture analyzers display to 0.001g or 0.01% moisture. When specifications call for very low moisture content such as for certain plastic resins select a balance with a resolution of 0.0001g or 0.001% moisture.
2. Repeatability - the ability of a moisture balance to display the same result when the same sample is tested repeatedly. It is expressed as a standard deviation.
3. Specialized capabilities such as providing graphic displays of the drying profile, and building libraries of frequently performed test parameters.
4. Temperature range – range from 50° to 160°C to broader ranges such as 30° to 200°C. The temperature can usually be set in increments of 1°C.
5. Some models have only one drying mode and others have several, such as standard, soft, fast and step.
6. Switch-off criteria by time, weight loss per unit of time or manual.
7. Maximum sample weight with 50 to 70g typical but some units have higher capacities.
8. Some models come with a calibration weight and for others the weight must be purchased separately.
9. Warranties vary.

### **How do I Decide on Optional Equipment?**

“Standard” and “optional” equipment are defined by the moisture analyzer manufacturer. Certain items described above such as temperature calibration sets are usually optional. The same applies to this list of accessories or options that will prove useful immediately or down the line:

- Aluminum sample plates
- Fiberglass filters
- Printers – matrix or thermal
- Calibration weights if not included
- RS 232 Interface if not included

We hope the information here is useful when selecting a moisture analyzer. For additional information and recommendations on selecting equipment please contact the scientists at Tovatech for expert advice on the right unit for your applications.

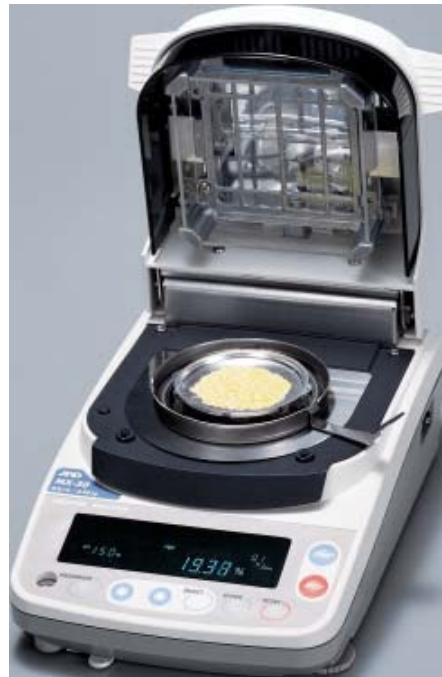

Need to talk to someone about moisture analysis?



**CALL US NOW TO GET YOUR ANSWER**  
**(973) 913-9734**



**CLICK HERE TO**  
**CHAT WITH US NOW**



**CLICK HERE TO**  
**WRITE TO US NOW**